↔ **DOMETIC**

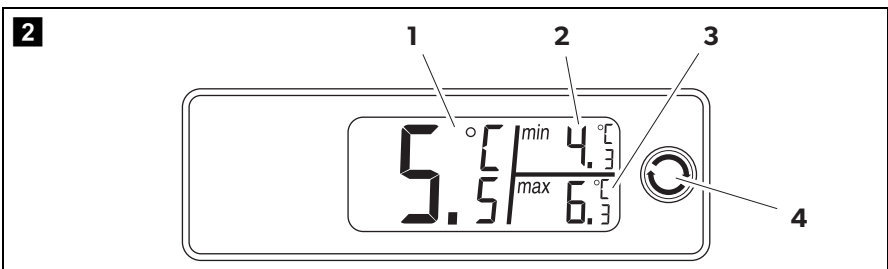
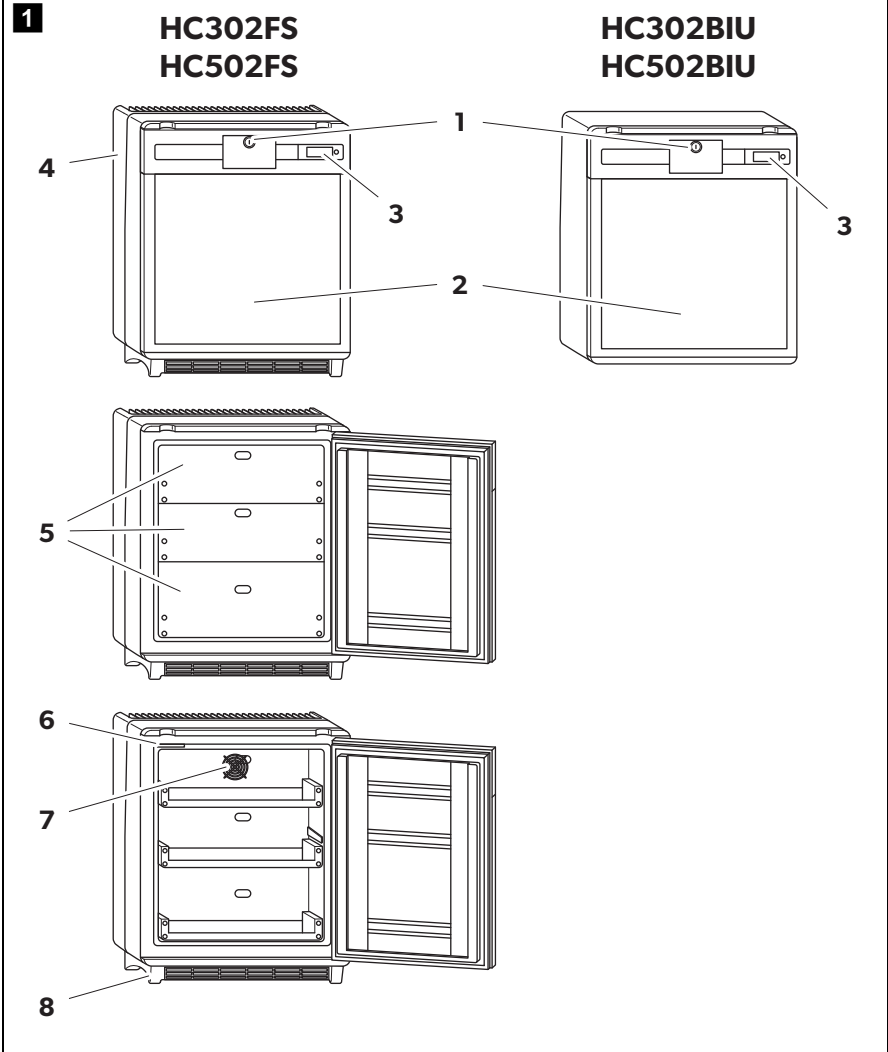
# PROFESSIONAL COOLING PHARMACEUTICAL

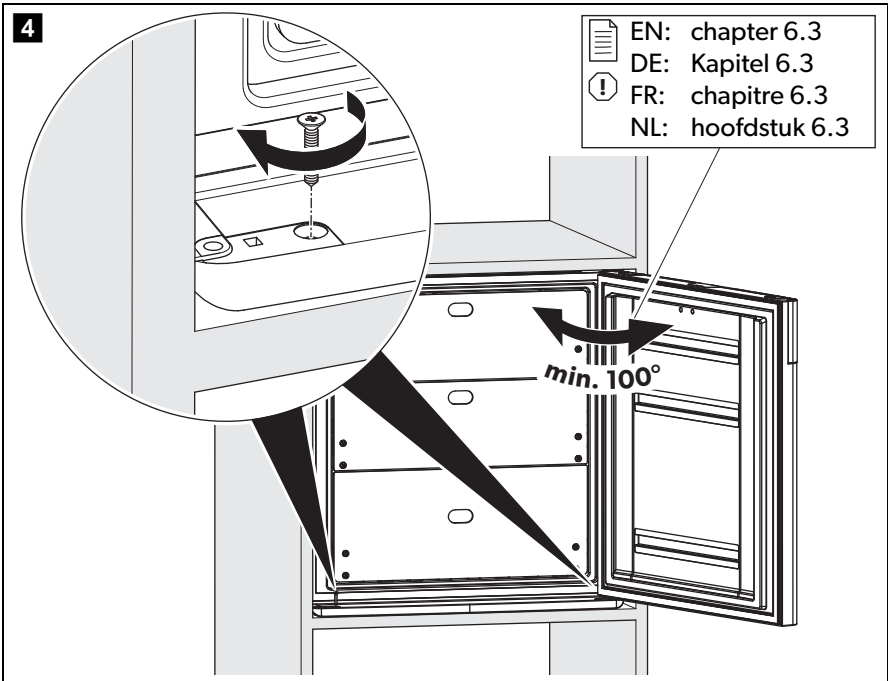
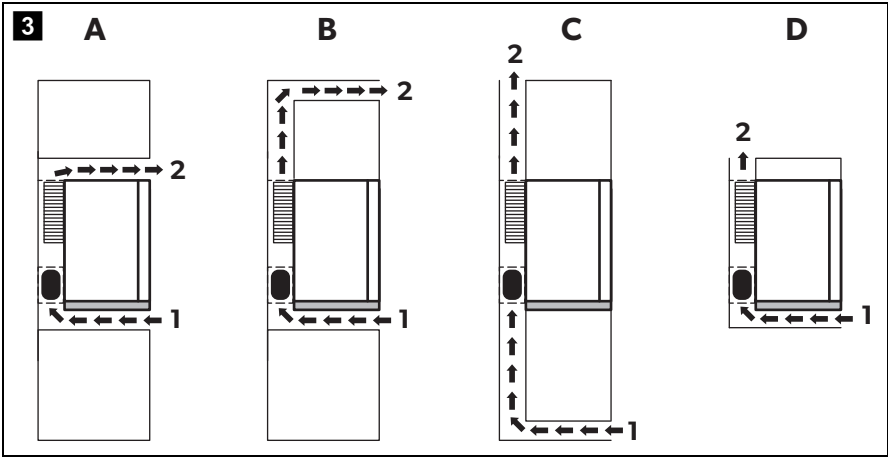
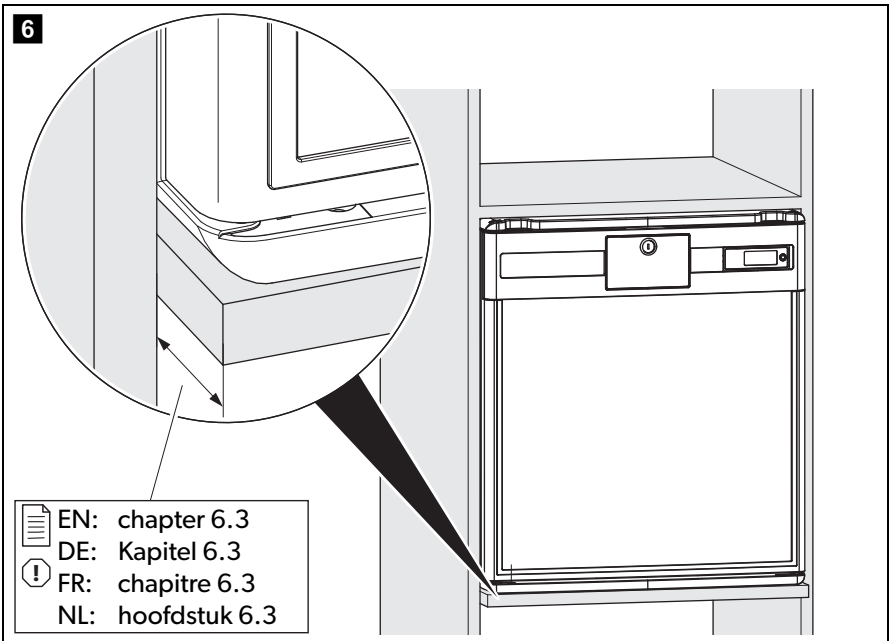
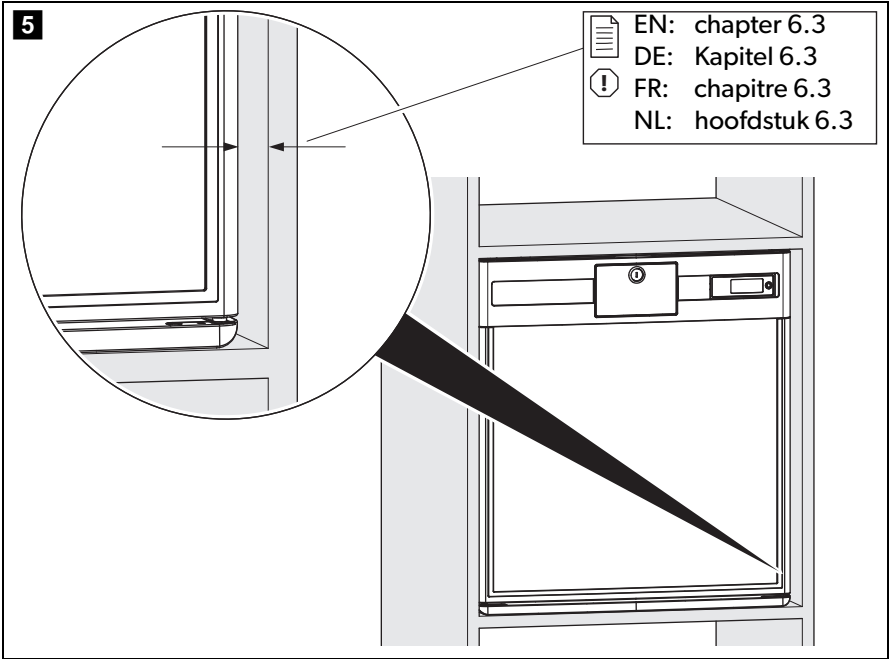
**HC302, HC502**

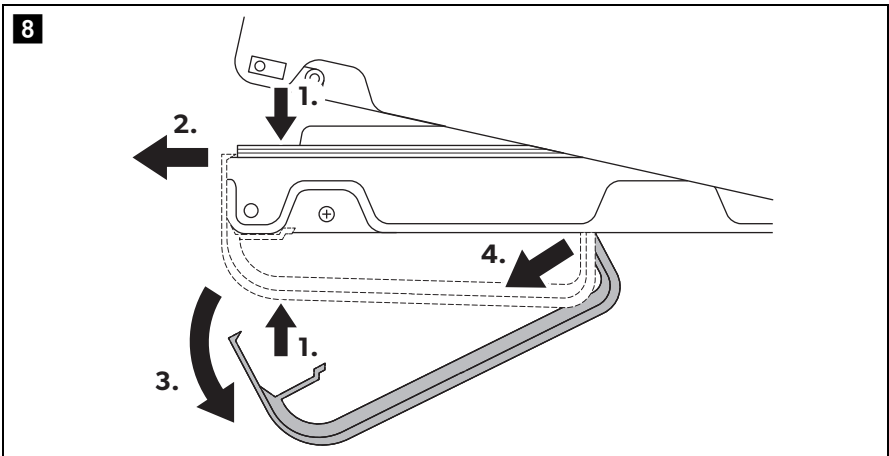
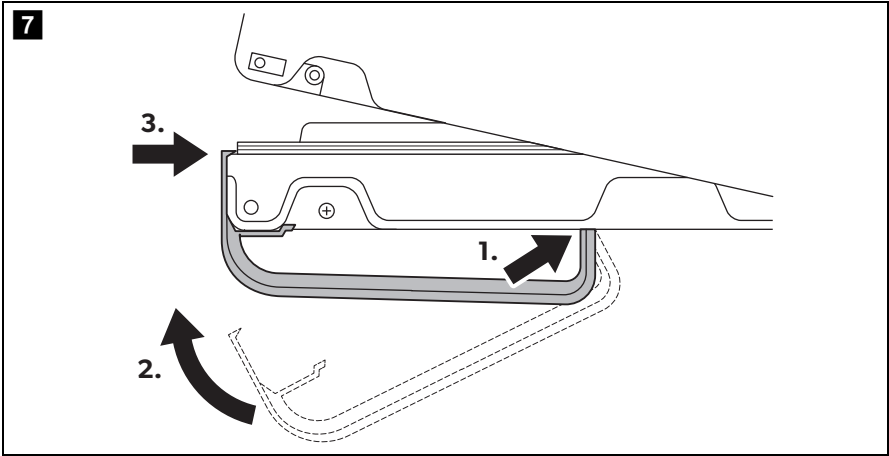
<b>EN</b>	<b>Health care refrigerator</b> Installation and Operating Manual . . . . .	9
<b>DE</b>	<b>Medikamentenkühlschrank</b> Montage- und Bedienungsanleitung . . . . .	23
<b>FR</b>	<b>Réfrigérateur pour médicaments</b> Instructions de montage et de service . . . . .	38
<b>NL</b>	<b>Medicamentenkoelkast</b> Montagehandleiding en gebruiksaanwijzing . . . . .	53

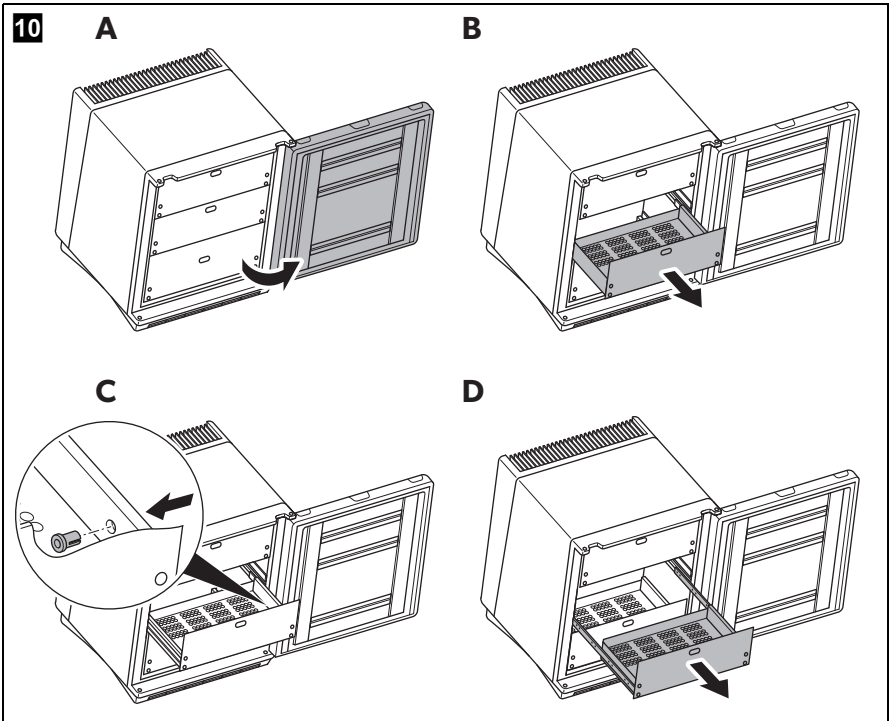
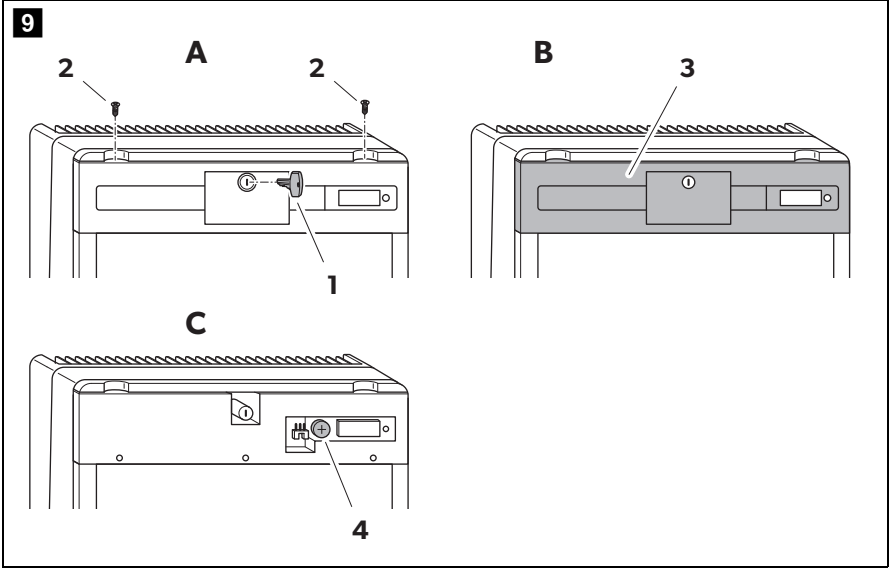
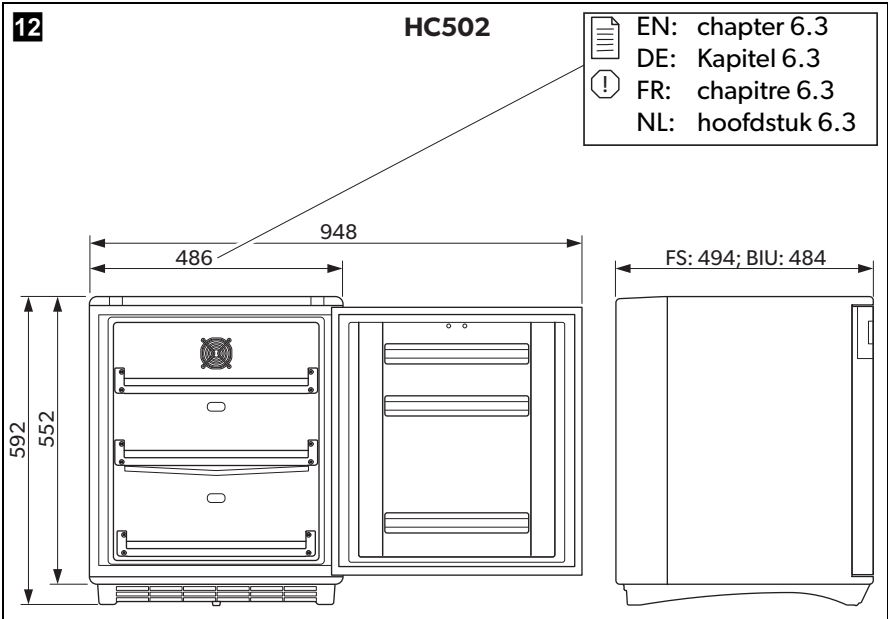
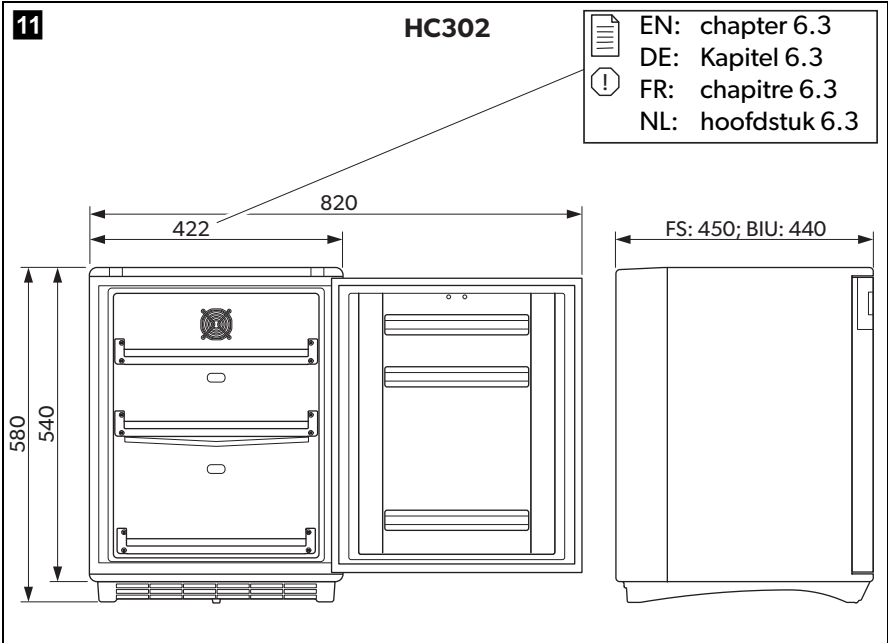
**Please read this instruction manual carefully before installation and first use, and store it in a safe place. If you pass on the product to another person, hand over this instruction manual along with it.**

## **Table of contents**

1	Explanation of symbols . . . . .	10
2	Safety instructions . . . . .	10
3	Scope of delivery . . . . .	12
4	Intended use . . . . .	13
5	Technical description . . . . .	13
6	Installing and connecting the refrigerator . . . . .	14
7	Using the refrigerator . . . . .	16
8	Cleaning and maintenance. . . . .	19
9	Troubleshooting . . . . .	20
10	Warranty . . . . .	21
11	Disposal . . . . .	21
12	Technical data . . . . .	22

# 1 Explanation of symbols

**DANGER!**

**Safety instruction:** Failure to observe this instruction will cause fatal or serious injury.

**WARNING!**

**Safety instruction:** Failure to observe this instruction can cause fatal or serious injury.

**CAUTION!**

**Safety instruction:** Failure to observe this instruction can lead to injury.

**NOTICE!**

Failure to observe this instruction can cause material damage and impair the function of the product.

**NOTE**

Supplementary information for operating the product.

## 2 Safety instructions

The manufacturer accepts no liability for damage in the following cases:

- Faulty assembly or connection
- Damage to the product resulting from mechanical influences and excess voltage
- Alterations to the product without express permission from the manufacturer
- Use for purposes other than those described in the operating manual

### 2.1 General safety

**WARNING!**

- Do not operate the device if it is visibly damaged.
- If this device's power cable is damaged, it must be replaced by customer service or a similarly qualified person in order to prevent safety hazards.
- This device should be repaired by qualified person only. Inadequate repairs may cause serious hazards.

- This device can be used by children aged 8 years or over, as well as by persons with diminished physical, sensory or mental capacities or a lack of experience and knowledge, providing they are supervised, or have been taught how to use the device safely and are aware of the resulting risks.
- Cleaning and user maintenance must not be carried out by children without supervision.
- Children must be supervised to ensure that they do not play with the device.
- Always keep and use the device out of the reach of children under 8 years of age.
- Do not store any explosive substances, such as spray cans with propellants, in the device.

**NOTICE!**

- Check that the voltage specification on the type plate is the same as that of the power supply.
- Never pull the plug out of the socket by the connection cable.
- The refrigerator is not suitable for storing substances which are caustic or contain solvents.
- Keep the drainage outlet clean at all times.
- Do not open the refrigerant circuit under any circumstances.
- Never transport the device in a horizontal position, so that no oil can leak out of the compressor.
- Make sure that the refrigerator circuit is not damaged during transportation.
- Set up the device in a dry location where it is protected against splashing water.

## 2.2 Operating the device safely

**DANGER!**

- Do not touch exposed cables.

**CAUTION!**

- Before starting the device, ensure that the power supply line and the plug are dry.

**NOTICE!**

- Do not use electrical devices inside the cooling device unless they are recommended by the manufacturer for that purpose.
- Do not place it near naked flames or other heat sources (heaters, direct sunlight, gas ovens etc.)
- **Danger of overheating!**  
Always ensure sufficient ventilation so that the heat generated during operation can dissipate. Make sure that the device is sufficiently far away from walls and other objects so that the air can circulate (fig. **3**, page 4).
- Ensure that the ventilation vents are not covered.
- The device shall not be exposed to rain.
- Never immerse the appliance in water.
- Protect the appliance and power cable against heat and moisture.
- Make sure that the stored substances do not touch the walls of the cooling area.

### 3 Scope of delivery

Quantity	Description
1	Refrigerator
1	Handle
1	Battery (CR1632)
1	Installation and operating manual

These instructions are also available in PDF format (e.g. on the manufacturer's website).

## 4 Intended use

The refrigerator is designed to store medicaments and medical products which may be stored at temperatures +2 °C and +8 °C.

**CAUTION!**

The refrigerator is not suitable for the storage of blood.

## 5 Technical description

### 5.1 Function description

All materials used in the refrigerator are compatible for use with medicines.

The refrigerant circuit is maintenance-free.

The refrigerator is available as freestanding version (FS) and as built-in version (BIU).

The freestanding version (HC 302FS, HC502FS) is intended for freestanding use. The cooling unit cover protects the warm parts of the cooling units. The foot frame with ventilation openings ensure sufficient ventilation. By removing the foot frame and cooling unit cover the refrigerator can be used as a built-in version.

The built-in version (HC 302BIU, HC502BIU) is intended to be mounted into furniture. It doesn't have a foot frame or cooling unit cover.

The digital thermometer shows and stores the minimum and maximum inner temperatures from the last reset so that you can check whether the medicines have been stored at the appropriate temperatures.

## 5.2 Control elements

No. in fig. <b>1</b> , page 3	Explanation
1	Door lock
2	Door
3	Digital thermometer (see fig. <b>2</b> , page 3)
4	Cooling unit cover
5	Drawers
6	Interior light
7	Fan
8	Foot frame

Key to fig. **2**, page 3

No.	Explanation
1	Current inner temperature (in °C)
2	Minimum inner temperature (in °C)
3	Maximum inner temperature (in °C)
4	Reset button

# 6 Installing and connecting the refrigerator

## 6.1 Unpacking the refrigerator

- Check that the packaging is not damaged.
- Remove the cardboard packaging.
- Check that the refrigerator is not damaged.

If you find damage on the packaging or at the refrigerator resulting from transport, report it immediately to the transportation firm.

## 6.2 Installing freestanding version

Observe the following instructions when installing:

- Install the refrigerator so that the warm air produced can easily flow away, either upwards or to the sides.
  - Leave the device to stand for at least four hours before you connect it to the power supply so that the refrigerant can flow back into the compressor.
- Place the refrigerator at the chosen place.

## 6.3 Installing built-in version

Observe the following instructions when installing:

- Install the refrigerator so that the warm air produced can easily flow away (either upwards or to the front, fig. **3**, page 4).  
Key to fig. **3**, page 4

No.	Explanation
1	Cold intake air
2	Hot waste air

- Leave the device to stand for at least four hours before you connect it to the power supply so that the refrigerant can flow back into the compressor.
  - Make sure that the refrigerator door is able to open at least 100°. Otherwise the drawers cannot be used.
- Mount the refrigerator to the furniture by two screws through the fittings at the bottom (fig. **4**, page 4).

## 6.4 Connecting the refrigerator to the power supply

### NOTICE!

**UK only:** The refrigerator is supplied with a plug fitted with 3 A fuse. In case this fuse has to be changed, a 3 A ASTA approved (BS 1362) fuse must be used.

- Connect the plug of the connection cable into the AC mains.
- Insert the battery into the thermometer (see chapter “Replacing the battery of the thermometer” on page 17).
- ✓ The display turns on.  
For the first two seconds the display will not show temperature values; it shows “--.”. Then the temperature read from the sensor is shown.

## 6.5 Mounting and removing the door handle

- Mount the handle as shown (fig. **7**, page 6).
- Remove the handle as shown (fig. **8**, page 6).

# 7 Using the refrigerator

### NOTE

- Before starting your new refrigerator for the first time, you should clean it inside and outside with a damp cloth for hygienic reasons (please also refer to the chapter “Cleaning and maintenance” on page 19).
- Always leave some space between the medicaments in the cabinet to ensure efficient and even cooling. If the drawers are packed too tightly there may be unacceptable temperature values.

## 7.1 Energy saving tips

- Choose a well ventilated location which is protected from direct sunlight.
- Do not open the refrigerator more often than necessary.
- Do not leave the doors open for longer than necessary.
- Defrost your refrigerator as soon as a layer of ice forms.

## 7.2 Using the refrigerator

### NOTE

The temperature values are shown in °C. This cannot be changed.

After the first start, after replacing the battery or after switching off the refrigerator for a longer time the display of the digital thermometer (fig. **2**, page 3) shows the current ambient temperature. When the inner temperature reaches the operating range the thermometer should be reset to avoid showing wrong maximum temperature value.

- Press the reset button (fig. **2** 4, page 3) for 5 s.

The thermometer has a blue backlight which makes the temperature values easier to read.

- Press the reset button (fig. **2** 4, page 3) briefly.
- ✓ The blue backlight lights for 2 s.

The thermometer stores the values of the minimum and maximum inner temperature since the last reset. These values are deleted when the thermometer is reset or the battery is changed.

## 7.3 Replacing the battery of the thermometer

Replace the battery of the thermometer as follows (fig. **9**, page 7):

- Pull the key (**1**) from the lock.
- Remove the lock cover rosette using two standard screwdrivers.
- Remove both screws (**2**) from front panel.
- Push the front panel (**3**) a bit upwards and remove it by slightly pulling.
- Remove the battery (**4**).

### NOTICE!

Make sure the polarity is correct (the “+” must face up).

- Insert the new battery (CR1632) in its holder.
- Remount the front panel and lock cover rosette in reversed order.

## 7.4 Defrosting the refrigerator

### NOTE

In high ambient temperatures, the refrigerator may operate continuously, which can lead to excessive formation of frost on the rear wall of the refrigerator compartment.

Humidity can cause frost to form in the interior of the refrigerator or on the vaporiser.

A slight layer of frost or water droplets could form in the interior of the refrigerator if it has been cooling for longer periods. This is normal because the moisture in the air condenses to water when the temperature in the refrigerator drops. The refrigerator is **not** defective.

The condensate flows out through a water drain duct into a water container at the back of the unit at top of the compressor, where it evaporates.

It is recommended to defrost the refrigerator every three weeks.

Proceed as follows:

- Disconnect the refrigerator from the AC mains.
- Remove the drawers (see chapter “Removing the drawers” on page 19).
- Leave the door open.
- After defrosting wiped out the cabinet interior with a clean, damp cloth.
- Assemble the drawers.

## 7.5 Switching off and storing the refrigerator

If you do not intend to use the refrigerator for a long time, proceed as follows:

- Disconnect the plug.
- Clean the refrigerator (see chapter “Cleaning and maintenance” on page 19).
- Leave the door slightly open.  
This prevents smells from building up.

## 7.6 Removing the drawers

You can remove the drawers, e.g for cleaning.

Proceed as follows (fig. **10**, page 7):

- Open the door (**A**).
- Pull out the drawer (**B**).
- Push the fixing pins from the outside to the inside (**C**)
- Remove the drawer (**D**)
- Assemble the drawers in reverse order.

## 8 Cleaning and maintenance

### WARNING!

Always disconnect the refrigerator from the mains before you clean and service it.

### NOTICE!

- Do not use abrasive cleaning agents or hard objects during cleaning as these can damage the refrigerator.
- Never use hard or sharp tools to remove ice or to free objects frozen onto the device.
- Do not use any mechanical tools or any other tools to speed up the defrosting process.

- Clean the refrigerator regularly and as soon as it becomes dirty with a damp cloth.
- Make sure that no water drips into the seals. This can damage the electronics.
- Wipe the refrigerator dry with a cloth after cleaning.
- Check the condensate drain regularly.  
Clean the condensate drain when necessary. If it is blocked the condensate collects on the bottom of the refrigerator.
- Regularly clean the door seals.
- **Freestanding version:** Clean the grid of the condenser and the compressor on the back of the device with a vacuum cleaner regularly.

# 9 Troubleshooting

## Water in or under the refrigerator

Fault	Possible cause	Remedy
Water is leaking into the inside of the refrigerator.	The condensate drain duct is clogged.	Clean the drain.
	Medicines are preventing the water from flowing to the collector.	Make sure that medicines do not touch the rear wall.
Water is leaking onto the floor.	There is too much water in the water container.	Clean the water container.

## Display signals

Signal	Explanation
<p><b>General error</b></p> <p>The display blinks each second.</p>	The display shows alternately the current inner temperature and the active error code (fig. 2 1, page 3) (E1 or E2, see below).
<p><b>Sensor failure</b></p> <p>The display blinks and shows "E1".</p>	The minimum and maximum values will be kept as the last known values.
<p><b>Low battery</b></p> <p>The display blinks and shows "E2" alternately with the current temperature.</p>	The minimum and maximum inner temperatures are still visible and will be updated from the current inner temperature reading.

## 10 Warranty

The statutory warranty period applies. If the product is defective, please contact the manufacturer's branch in your country (see the back of the instruction manual for the addresses) or your retailer.

For repair and guarantee processing, please include the following documents when you send in the device:

- A copy of the receipt with purchasing date
- A reason for the claim or description of the fault

## 11 Disposal

### **WARNING! Children beware!**

Before disposing of your old refrigerator:

- Dismantle the drawers.
- Leave the drawers in the refrigerator so that children cannot climb inside.

### **NOTICE!**

An insulating gas is used in the device (see type plate). Check with the relevant local government authority or your local waste disposal company about proper disposal and make sure that the device and all materials are disposed of professionally.

- Place the packaging material in the appropriate recycling waste bins wherever possible.

If you wish to finally dispose of the product, ask your local recycling centre or specialist dealer for details about how to do this in accordance with the applicable disposal regulations.

## 12 Technical data

	HC 302	HC 502
Gross capacity:	33 l	49 l
Capacity:	29 l	43 l
Connection voltage:	220 – 240 V	
Energy consumption:	0.42 kWh/24 h	0.44 kWh/24 h
Power consumption:	0.54 A	
Cooling capacity:	+2 °C to +8 °C	
Climatic class:	SN	
Refrigerant:	R134a	
Refrigerant quantity:	38 g	
CO <sub>2</sub> equivalent:	0.054 t	
Global warming potential (GWP):	1430	
Dimensions:	see fig. <b>11</b> , page 8	see fig. <b>12</b> , page 8
Weight:	approx. 20 kg	approx. 22 kg
Inspection/certification:		

Complies with the following EEC directives:

- LVD directive: 2006/95/EC
- EMC directive: 2004/108/EC
- CE directive: 93/68/EC
- RoHS directive: 2002/95/EC
- WEEE directive: 2002/96/EC

Contains fluorinated greenhouse gases

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**GERMANY****Dometic WAECO International GmbH**

Hollefeldstraße 63 · D-48282 Emsdetten  
 ☎ +49 (0) 2572 879-195 · 📠 +49 (0) 2572 879-322  
 Mail: info@dometic-waeco.de

**dometic.com****AUSTRALIA****Dometic Australia Pty. Ltd.**

1 John Duncan Court  
 Varsity Lakes QLD 4227  
 ☎ 1800 212121  
 📠 +61 7 55076001  
 Mail: sales@dometic.com.au

**AUSTRIA****Dometic Austria GmbH**

Neudorferstraße 108  
 A-2353 Guntramsdorf  
 ☎ +43 2236 908070  
 📠 +43 2236 90807060  
 Mail: info@dometic.at

**BENELUX****Dometic Branch Office Belgium**

Zincstraat 3  
 B-1500 Halle  
 ☎ +32 2 3598040  
 📠 +32 2 3598050  
 Mail: info@dometic.be

**BRAZIL****Dometic DO Brasil LTDA**

Avenida Paulista 1754, conj. 111  
 SP 01310-920 Sao Paulo  
 ☎ +55 11 3251 3352  
 📠 +55 11 3251 3362  
 Mail: info@dometic.com.br

**DENMARK****Dometic Denmark A/S**

Nordensvej 15, Taulov  
 DK-7000 Fredericia  
 ☎ +45 75585966  
 📠 +45 75586307  
 Mail: info@dometic.dk

**FINLAND****Dometic Finland OY**

Mestariitie 4  
 FIN-01730 Vantaa  
 ☎ +358 20 7413220  
 📠 +358 9 7593700  
 Mail: info@dometic.fi

**FRANCE****Dometic SAS**

ZA du Pré de la Dame Jeanne  
 B.P. 5  
 F-60128 Pleilly  
 ☎ +33 3 44633525  
 📠 +33 3 44633518  
 Mail: vehiculesdeloisirs@dometic.fr

**HONG KONG****Dometic Group Asia Pacific**

Suites 2207-11 / 22/F - Tower 1  
 The Gateway - 25 Canton Road,  
 Tsim Sha Tsui - Kowloon  
 ☎ +852 2 4611386  
 📠 +852 2 4665553  
 Mail: info@waeco.com.hk

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**NEW ZEALAND****Dometic New Zealand Ltd.**

PO Box 12011  
 Penrose  
 Auckland 1642  
 ☎ +64 9 622 1490  
 📠 +64 9 622 1573  
 Mail: customerservices@dometic.co.nz

**NORWAY****Dometic Norway AS**

Østerøyveien 46  
 N-3232 Sandefjord  
 ☎ +47 33428450  
 📠 +47 33428459  
 Mail: firmapost@dometic.no

**POLAND****Dometic Poland Sp. z o.o.**

Ul. Puławska 435A  
 PL-02-801 Warszawa  
 ☎ +48 22 414 3200  
 📠 +48 22 414 3201  
 Mail: info@dometic.pl

**PORTUGAL****Dometic Spain, S.L.**

Branch Office em Portugal  
 Rot. de São Gonçalo nº 1 – Esc. 12  
 2775-399 Carcavelos  
 ☎ +351 219 244 173  
 📠 +351 219 243 206  
 Mail: info@dometic.pt

**RUSSIA****Dometic RUS LLC**

Komsomolskaya square 6-1  
 RU-107140 Moscow  
 ☎ +7 495 780 79 39  
 📠 +7 495 916 56 53  
 Mail: info@dometic.ru

**SINGAPORE****Dometic Pte Ltd**

18 Boon Lay Way 06-140 Trade Hub 21  
 Singapore 609966  
 ☎ +65 6795 3177  
 📠 +65 6862 6620  
 Mail: dometic@dometic.com.sg

**SLOVAKIA****Dometic Slovakia s.r.o. Sales Office Bratislava**

Nádražná 34/A  
 900 28 Ivanka pri Dunaji  
 ☎/📠 +421 2 45 529 680  
 Mail: bratislava@dometic.com

**SOUTH AFRICA****Dometic (Pty) Ltd.****Regional Office****South Africa & Sub-Saharan Africa**

2 Avalon Road  
 West Lake View Ext 11  
 Modderfontein 1645  
 Johannesburg  
 ☎ +27 11 4504978  
 📠 +27 11 4504976  
 Mail: info@dometic.co.za

**SPAIN****Dometic Spain S.L.**

Avda. Sierra del Guadarrama, 16  
 E-28691 Villanueva de la Cañada  
 Madrid  
 ☎ +34 91 833 60 89  
 📠 +34 900 100 245  
 Mail: info@dometic.es

**SWEDEN****Dometic Scandinavia AB**

Gustaf Melins gata 7  
 S-42131 Västra Frölunda  
 ☎ +46 31 7341100  
 📠 +46 31 7341101  
 Mail: info@dometicgroup.se

**SWITZERLAND****Dometic Switzerland AG**

Riedackerstrasse 7a  
 CH-8153 Rümlang  
 ☎ +41 44 8187171  
 📠 +41 44 8187191  
 Mail: info@dometic.ch

**UNITED ARAB EMIRATES****Dometic Middle East FZCO**

P. O. Box 17860  
 S-D 6, Jebel Ali Freezone  
 Dubai  
 ☎ +971 4 883 3858  
 📠 +971 4 883 3868  
 Mail: info@dometic.ae

**UNITED KINGDOM****Dometic UK Ltd.**

Dometic House, The Brewery  
 Blandford St. Mary  
 Dorset DT11 9LS  
 ☎ +44 344 626 0133  
 📠 +44 344 626 0143  
 Mail: customerservices@dometic.co.uk

**USA****Dometic RV Division**

1120 North Main Street  
 Elkhart, IN 46515  
 ☎ +1 574-264-2131