Istruzioni per l'utilizzatore
User Instruction
Instructions pour l'utilisateur
Benutzeranleitungen

ARMADIO REFRIGERATO
REFRIGERATED CABINET
ARMOIRE REFRIGEREE
KÜHLSCHRANK

xPRO

VERS. 0816

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ENGLISH**Content****8 USER INSTRUCTIONS**

8.1 CONTROLS	17
8.1.1 DESCRIPTION OF CONTROLS AND TOUCH-BUTTONS OF THE xPRO CONTROL PANEL	17
8.1.2 THE DISPLAY AND THE FUNCTIONING INDICATIONS	17
8.1.3 AUTOMATIC/MANUAL TOUCH-BUTTONS SAFETY LOCKING PROTECTION AND UNLOCKING CODE	18
8.2 OPERATION OF THE xPRO CONTROL PANEL	18
8.2.1 START-UP.....	18
8.2.2 SEQUENCE AND ON/OFF CODE	18
8.2.3 STOPPING THE UNIT	18
8.2.4 DATE AND TIME SETTING	19
8.2.5 GENERAL INFORMATION OF THE xPRO CONTROL PANEL FUNCTIONING	19
8.3 THE MAIN MENU	19
8.3.1 SETPOINT MENU	19
8.3.2 DATAS MENU	20
8.3.2.1 Temperature chart	20
8.3.2.2 Battery level	20
8.3.2.3 Functioning timers	20
8.3.2.4 Door openings	20
8.3.2.5 Defrosting	21
8.3.2.6 Sensors	21
8.3.3 MANUAL DEFROST	21
8.3.4 ALARMS MEMORY	21
8.3.5 GENERAL TEST AND TEST FOR TEMPERATURES ALARM	21
8.3.6 SETTING MENU	22
8.3.6.1 Language setting	22
8.3.6.2 Alarm setting	23
8.3.6.3 Ventilation setting	24
8.3.6.4 Download USB data	24
8.3.6.5 Reactivation alarm after muting	24
8.3.6.6 Temperature alarm activation delay	24
8.3.6.7 Door ajar alarm activation delay	25
8.3.6.8 Codes setting	25
8.3.6.9 Calibration	25
8.4 OPERATION OF THE TEMPERATURE CHART RECORDER	26
8.4.1 DESCRIPTION OF THE TEMPERATURE CHART RECORDER	26
8.4.1.1 Paper chart replacement	26
8.4.1.2 Ink-tip replacement	26
8.4.1.3 Battery replacement	26
8.5 OPERATION OF THE DIGITAL PRINTER	27
8.5.1 DESCRIPTION OF THE DIGITAL PRINTER	27

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8 USER INSTRUCTIONS

The information in this section of the manual regards the user or other non-specialized personnel (see par. 1.3 in the “Instruction and Maintenance Manual”).

After the appliance has been installed in accordance with the instructions of section 3 of the “Instruction and Maintenance Manual”, it is ready for use.

8.1 CONTROLS

According to the models, the appliance is equipped of different types of controls:

- xPRO control panel (*Fig.10*)
- Temperature chart recorder (*Fig.22*)

8.1.1 Description of controls and touch-buttons of the xPRO control panel (*Fig.10*)

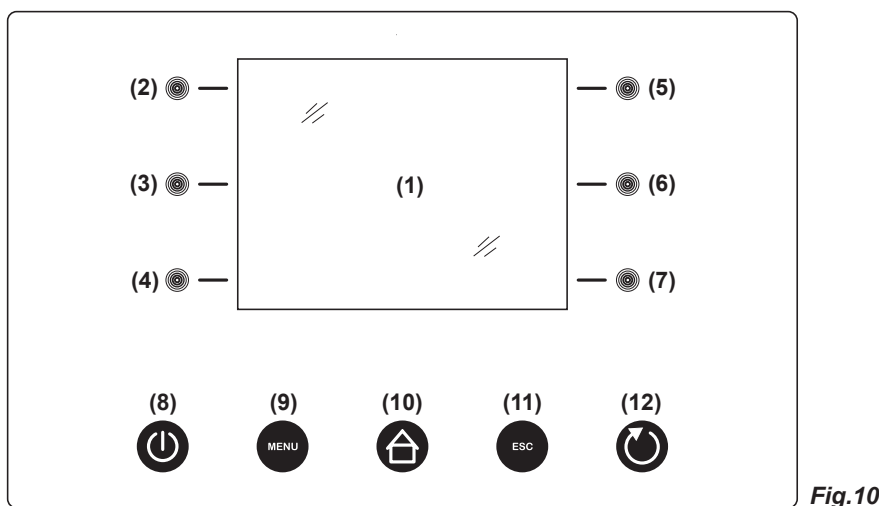
Fig.10

The appliance is equipped with the following controls:

- Display (1). Showing and controlling the temperature and the various machine functioning states
- Capacitive **touch-buttons** (2), (3), (4), (5), (6), (7). Allow various functions according to the OSD (on screen display) menu shown on the display
- **ON/OFF touch-button** (8). Allows the display/appliance switching ON and OFF
- **MENU touch-button** (9). Allows to enter into the main menu
- **HOME touch-button** (10). Allows (if back-lightened) to return back to the home (main) screen-shot
- **ESC touch-button** (11). Allows to return back to the previous menu
- **ENTER touch-button** (12). Allows to enter/confirm the variations made on the function/menu shown on the display

8.1.2 The display (1) and the functioning indications (*Fig.11*)

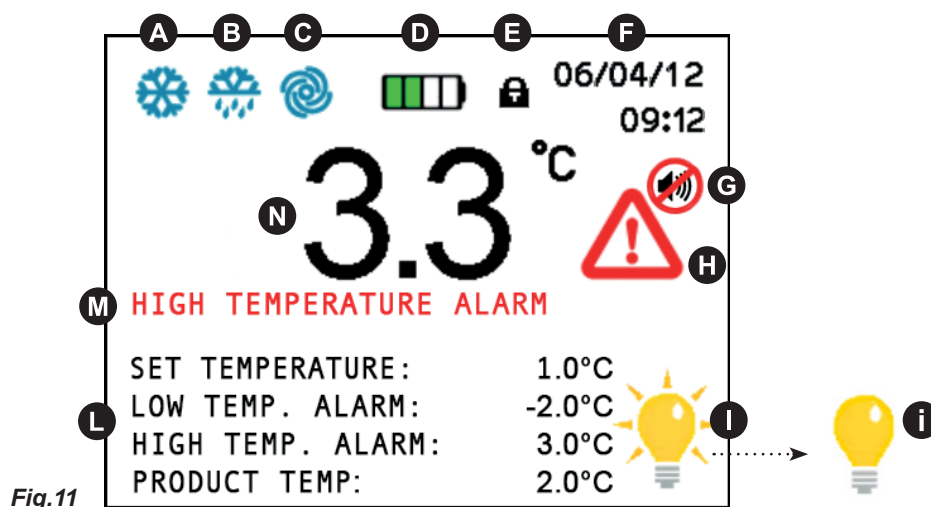
Fig.11

The display shows the below listed icons/indications that, if enlightened, give indications/alert of the running functions:

- **A** COMPRESSOR: lights steadily when the compressor is running, blinks in case of starting delay or when protection is active
- **B** DEFROST: lights steadily during defrosting cycle

- **C** FAN: lights steadily when the evaporator fan is working
- **D** BATTERY: shows the actual charging status of the back-up battery or if a e-charging is in progress
- **E** TOUCH-BUTTONS LOCKING: appears and lights steadily when the touch-buttons are in protection locking mode
- **F** DATE/TIME: shows the actual date and time
- **G** BUZZER: lights steadily when acoustic alarm indication has been muted through the **touch-button (6)**
- **H** ALARM ALERT: lights steadily when an alarm is in progress. This icon is always accompanied from the relative alarm description **M**
- **I** INTERNAL LIGHTING: lights steadily when the internal lighting is activated. Pressing the **touch-button (7)**, it is possible to switch manually the internal lighting ON or OFF. When the internal lighting is not activated, it appears the icon **i**. The internal lighting switch also ON automatically at every door opening and switch OFF at the closing.
- **L** SETTINGS: show the actual appliance settings, set temperature, high and low alarm limits and if present (optionally available) the actual temperature of the stored product (simulation)
- **N** ACTUAL TEMPERATURE: shows the actual internal temperature of the appliance. In case the temperature is out of the allowed limits, the indication of the temperature value pass to red colour

8.1.3 Automatic/manual touch-buttons safety locking protection and unlocking code

The xPRO control panel touch-buttons are provided with a safety locking protection to avoid tampering from unauthorized personnel. The key locking protection has an automatic activation after 5 minutes from the last operation executed with the touch-buttons. The key locking condition is signalled on the display, top part, with the icon of the lock-pad **E**. To re-enable the use of the touch-buttons, press any touch-button and enter the **unlocking code 0010**. To activate manually the touch-buttons safety locking protection, press the **ESC touch-button (11)** and the **HOME touch-button (10)** even if not enlightened.

8.2 OPERATION OF THE xPRO CONTROL PANEL

8.2.1 Start-up

Before starting up the unit check that the electrical connections have been made correctly as indicated in headings 3.3 and 3.4 "Instruction and Maintenance Manual".

Perform preliminary cleaning of the unit as described in sub-heading 5.2.1 "Instruction and Maintenance Manual".

At the starting, when the xPRO control panel is powered, it will appear a screen-shot with the EVERmed logo for approx. 10 seconds. Immediately after, the display will show the STAND-BY condition with the date and time with white indications on a black back-ground.

8.2.2 Sequence and ON/OFF code (Fig.10)

- if the appliance is switched OFF, but connected to the electrical network, the display shows the STAND-BY condition with the date and time with white indications on a black back-ground
- press the **ON/OFF touch-button (8)** to switch the appliance ON; the display shows the screen-shot to insert the ON/OFF code
- enter the **ON/OFF code 0050** through the **touch-buttons (3) (6) and (4) (7)**
- the touch-buttons **(3)←** and **(6)→** have to be used to select the position where to enter the numerical value indicated from the cursor that can be moved right and left
- the touch-buttons **(4)-** and **(7)+** have to be used to lower and/or increase the highlighted numerical value
- having inserted the correct code, press the **ENTER touch-button (12)** to confirm, the appliance switch-on and the access to all touch-buttons is enabled

At every appliance starting, it will appear the indication "Battery level low" which has to be reset pressing the enlightened touch-button **(6)**. The battery low level alarm that appear at the starting will not be memorized in the alarms memory as it is not a real alarm, but it is only the indication of the battery-charger starting.

8.2.3 Stopping the unit (Fig.10)

- deactivate the eventual keyboard touch-buttons locking protection (see par. 8.1.3)
- with control panel on, press the **ON/OFF touch-button (8)** to switch the appliance OFF
- the display shows the screen-shot to insert the ON/OFF code
- enter the **ON/OFF code 0050 through the touch-buttons (3) (6) and (4) (7)**
- the touch-buttons **(3)←** and **(6)→** have to be used to select the position where to enter the numerical value indicated from the cursor that can be moved right and left
- the touch-buttons **(4)-** and **(7)+** have to be used to lower and/or increase the highlighted numerical value
- having inserted the correct code, press the **ENTER touch-button (12)** to confirm, the appliance switch-off

For further information about codes, see the menu CODE SETTING (par. 8.3.6.8).

8.2.4 Date and time setting (Fig.10)

To set the calendar and clock proceed as follow:

- deactivate the eventual touch-buttons locking protection (see par. 8.1.3)
- press the **MENU touch-button (9)** to access to the main menu (Fig.12)
- press the **touch-button (7)** to access to the setting menu (Fig.15)
- press the **touch-button (2)** to select the menu "Date and time"
- using the **touch-buttons (3)←** and **(6)→** highlight the position in the calendar or in the clock that should be modified
- using the **touch-buttons (4)-** and **(7)+** to set the desired value
- repeat the last 2 operations for all the numerical values in both calendar and clock, then press the **ENTER touch-button (12)** to confirm the new settings

8.2.5 General information of the xPRO control panel functioning

The xPRO control panel surveys and has the control of all the appliance functioning parameters and devices. The PRO control panel is provided with a back-up battery that grants its operations even in case of power failure, for about 48 (depending from the charge status). The battery charge status is automatically controlled through a special battery charging module in order to grant constantly the maximum battery efficiency. In the top part of the main screen-shot, it is always present an icon **D** indicating the battery status (showing charging in progress or the actual status of charge through the number of green colour charging steps).

The xPRO control panel alerts in case of abnormal functioning status with a both a visual indication (red colour icon **H**) and alarm description **M**) and an acoustic signalling through a buzzer.

The acoustic alarm has muting facilities, pressing the **touch-button (6)** (with icon **G** on display) and it is provided of automatic re-activation facilities once passed the pre-set time programmable through the specific menu "Reactivation alarm after muting" in case the alarm condition persists.

For further information about the alarms and their functioning check the paragraphs describing every single signalled alarm condition. Par. 8.3.4 and Par. 8.3.5.

8.3 THE MAIN MENU

To access to the Main Menu (Fig.12) press the **MENU touch-button (9)** (Fig.10). (enter the touch-buttons unlocking protection code **0010** if needed)

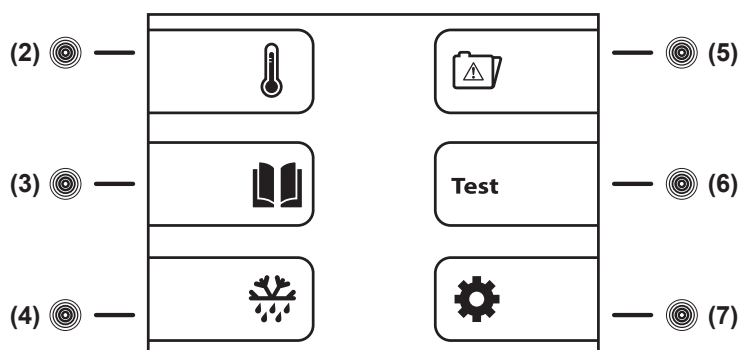
Fig.12

8.3.1 Setpoint Menu (touch-button (2) Fig.12)

The appliance is factory set for operation at the following temperature:

- serie MPR xPRO: +2°C / +15°C
- serie LR xPRO: 0°C / +15°C
- serie BBR xPRO: +4C
- serie LFG xPRO: -5°C / -20°C
- serie LF xPRO: -5°C / -20°C ~ -25°C
- serie LDF xPRO: -15°C / -30°C
- serie PDF xPRO: -25°C / -40°C
- serie LCRR xPRO: 0°C / +15°C (refrigerator compartment), 0°C / +15° (refrigerator compartment)
- serie MPRR xPRO: 2°C / +15°C (refrigerator compartment), 2°C / +15°C (refrigerator compartment)
- serie LCRF xPRO: 0°C / +15°C (refrigerator compartment), -5°C / -20°C ~ -25°C (freezer compartment)

To set the desired working temperature, proceed as follow:

- from the main menu screen-shot, press the **touch-button (2)** to access to the SETPOINT menu
- using the **touch-buttons (4)-** and **(7)+** to set the desired value
- press the **ENTER touch-button (12)** to confirm the new setting and the **ESC touch-button (11)** to return back to the previous menu

Press the **MENU touch-button** to return back to the main menu screen-shot without making any modification to the Setpoint value. If no touch-buttons are pressed for 50 seconds, automatically, the xPRO control panel return to the main screen-shot.

8.3.2 Datas Menu (touch-button (3) Fig.12)

From the main menu screen-shot (Fig.12), press the **touch-button (3)** to access to the Datas Menu. The Datas Menu appears as follow (Fig.13)

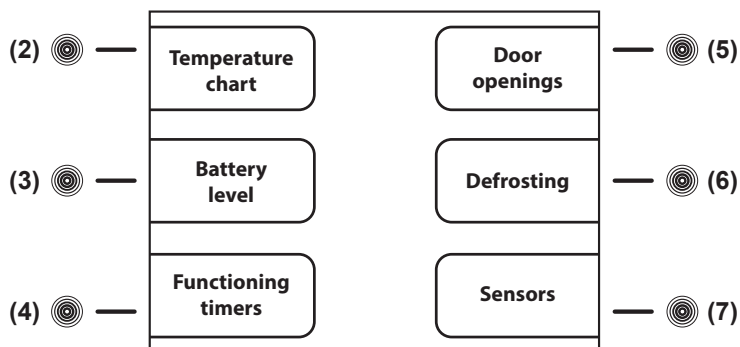
Fig.13

8.3.2.1 Temperature chart (touch-button (2) Fig.13)

The chart shows, in a graphical form, the temperature trend of the last operating period of the appliance.

- from the datas menu, press the **touch-button (2)** to access to the TEMPERATURE CHART menu

It is possible to operate on the chart displaying using the following touch-buttons:

- press the **touch-button (2)Y+** and the **touch-button (4)Y-**, the chart moves on the vertical ordinate axis, the temperature range, granting a wider view of the temperatures
- press the **touch-button (5)X+** and the **touch-button (7)X-**, the chart moves on the horizontal x-axis, allowing a more detailed monitoring of the chart trend in the time (till 600 minutes) and then in hours (till 46 hours)
- press the **touch-button (3)IN** and the **touch-button (6)OUT**, the chart zoom in and out for a enlarged or dwarfed view of the chart in its totality
- press the **ESC touch-button (11)** to return back to the previous menu or **HOME touch-button (10)** to return back to the main screen-shot

8.3.2.2 Battery level (touch-button (3) Fig.13)

To display the actual status of charge of the back-up battery, proceed as follow:

- from the datas menu, press the **touch-button (3)** to access to the BATTERY LEVEL screen-shot
- the display will show the word "CHARGING" if the battery is being charged or the number "100%" if the battery is fully charged
- press the **ESC touch-button (11)** to return back to the previous menu or **HOME touch-button (10)** to return back to the main screen-shot

8.3.2.3 Functioning timers (touch-button (4) Fig.13)

The functioning timers show important information about the appliance life and its working. This details can not be erased.

- from the datas menu, press the **touch-button (4)** to access to the FUNCTIONING TIMERS menu
- press the **touch-button (5)↑** and the **touch-button (7)↓**, it is possible to display the following timers:
 - Time Comp ON (average functioning times with compressor ON)
 - Time Comp OFF (average times with compressor OFF)
 - Time Comp2 ON (average functioning times with back-up compressor ON - if present)
 - Time Comp2 OFF (average times with back-up compressor OFF - if present)
 - % Comp ON per day (daily percentage of compressor functioning)
 - % Comp ON per hour (hourly percentage of compressor functioning)
 - % Comp2 ON per day (daily percentage of back-up compressor functioning - if present)
 - % Comp2 ON per hour (hourly percentage of back-up compressor functioning - if present)
 - Total hours Comp ON (total hours of compressor functioning)
 - Total hours Comp2 ON (total hours of back-up compressor functioning - if present)
 - Total hours appliance (appliance life)
- press the **ESC touch-button (11)** to return back to the previous menu or **HOME touch-button (10)** to return back to the main screen-shot

8.3.2.4 Door openings (touch-button (5) Fig.13)

It is possible to check the door events of the latest 50 days of the appliance functioning. For each day of functioning of the appliance, it is dedicated a page in the door log in which are indicated the total number of the door openings, how many of this have given an alarm and the total time that the door has been left opened.

To check the door opening log, proceed as follow:

- from the datas menu, press the **touch-button (5)** to access to the DOOR OPENINGS menu
- press the **touch-button (5)↑** and the **touch-button (7)↓** to scroll the daily reports of the door openings
- press the **ESC touch-button (11)** to return back to the previous menu or **HOME touch-button (10)** to return back to the main screen-shot

8.3.2.5 Defrosting (touch-button (6) Fig.13)

All the appliance defrosting are recorded in the defrosting log; on the display are shown the following details:

- date and time of the defrost starting
- the duration of the defrost
- the type of defrost, which is indicated from one of the following codes:
 - 01 - defrost started manually from the specific menu
 - 02 - defrost started after 1 hour from the appliance switch-ON
 - 03 - defrost started after a pre-fix working time of the appliance
 - 04 - defrost started after a pre-fix working time of the compressor
 - 06 - defrost started from automatic defrost
 - 07 - defrost started because of low evaporator temperature
 - 09 - defrost started from an alarm
- progressive recording number of the defrost

To check the defrosting log, proceed as follow:

- from the datas menu, press the **touch-button (6)** to access to the DEFROSTING menu
- press the **touch-button (5)↑** and the **touch-button (7)↓**, to scroll the various defrost reports
- press the **ESC touch-button (11)** to return back to the previous menu or **HOME touch-button (10)** to return back to the main screen-shot

8.3.2.6 Sensors (touch-button (7) Fig.13)

The screen-shot sensors allows to read the actual values of the various sensors installed in the appliance (storage chamber, evaporator, condenser and product simulation if present).

To check the sensor readings, proceed as follow:

- from the datas menu, press the **touch-button (7)** to access to the SENSORS menu
- press the **ESC touch-button (11)** to return back to the previous menu or **HOME touch-button (10)** to return back to the main screen-shot

8.3.3 Manual defrost (touch-button (4) Fig.12)

To process a manual defrost, proceed as follow:

- from the main menu screen-shot, press the **touch-button (4)** to access to the MANUAL DEFROST menu
- press the **touch-button (4)YES**, to process the manual defrost with visual indication in the main screen-shot through the related icon **B**
- press the **touch-button (7)NO**, to exit without process the defrost and return back to the main screen-shot

8.3.4 Alarms memory (touch-button (5) Fig.12)

With this menu is possible to check the latest 30 alarm conditions (and related informations) occurred to the appliance.

To check the alarm log, proceed as follow:

- from the main menu screen-shot, press the **touch-button (5)** to access to the ALARMS MEMORY menu
- press the **touch-button (5)↑** and the **touch-button (7)↓** to scroll the alarm log
- press the **ESC touch-button (11)** to return back to the previous menu or **HOME touch-button (10)** to return back to the main screen-shot

NB: For the alarms “Door ajar”, “High temperature”, “Battery level low” and “Expansion Board” it will be recorded even the highest peak of temperature reached into the appliance during the alarm condition.

For the alarms “Low temperature” it will be recorded even the lowest peak of temperature reached into the appliance during the alarm condition.

If the alarm condition is still in progress, it appears “in progress”.

In case there are no alarms recorded, the display shows NO DATA.

8.3.5 General test and test for temperatures alarm (touch-button (6) Fig.12)

To process the testing, proceed as follow:

- from the main menu screen-shot, press the **touch-button (6)** to access to the TEST menu (**Fig.14**)

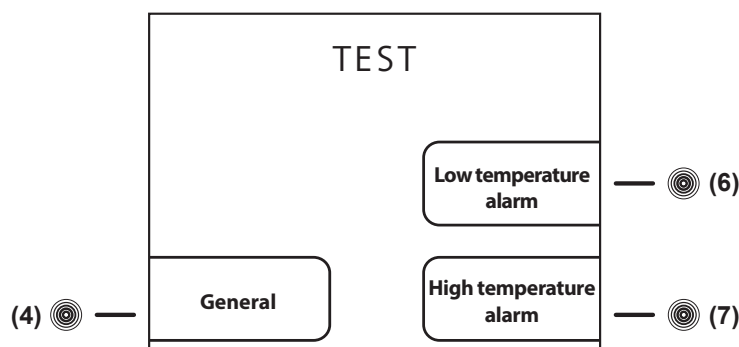
Fig.14

- press the **touch-button (4)** in the TEST screen-shot (**Fig.14**) to activate the GENERAL test. Through this procedure will be checked automatically the correct functioning of the battery and battery relay. The buzzer sounds during all the test (6 seconds). If the test has a positive result, the display shows TEST OK, otherwise the display switch off for a few seconds and automatically re-starts. Once completed the test, automatically, the display return back to the main screen-shot.

Once the appliance has reached the set temperature, it is possible to make the test temperature. Proceed as follows:

- press the **touch-button (6)** in the TEST screen-shot (**Fig.14**) to activate the LOW ALARM TEMPERATURE TEST. The display shows automatically the temperature that decrease and once reached the limit set for the low alarm temperature, It activates the buzzer and on the display appears the indication of Low temperature alarm. Press the **ESC touch-button (11)**, the test terminates and the display return back to the main screen-shot.

- press the **touch-button (7)** in the TEST screen-shot (**Fig.14**) to activate the HIGH ALARM TEMPERATURE TEST. The display shows automatically the temperature that increase and once reached the limit set for the high alarm temperature, It activates the buzzer and on the display appears the indication of Low temperature alarm. Press the **ESC touch-button (11)**, the test terminates and the display return back to the main screen-shot.

NB: During the test, the appliance continue to work normally and the alarm indication is not recorded in the alarms memory.

NB: In case of power blackout, the power is turned by the internal battery. The display will show the red message "**POWER FAILURE ALARM**" and the machine automatically shifts in the energy conservation mode. This status is recognized by the intermittently lighting of the display. To silence the buzzer momentarily press the touch (6). To silence the buzzer finally set the value "0" in the "Reactivation after muting" (Paragraph 8.3.6.5). This procedure will disable the buzzer reactivation of any subsequent alarms.

8.3.6 Setting menu  (touch-button (7) Fig.12)

To access to the SETTING menu, press the **touch-button (7)** from the main menu screen-shot. The access is to the first page of the menu (**Fig.15**).

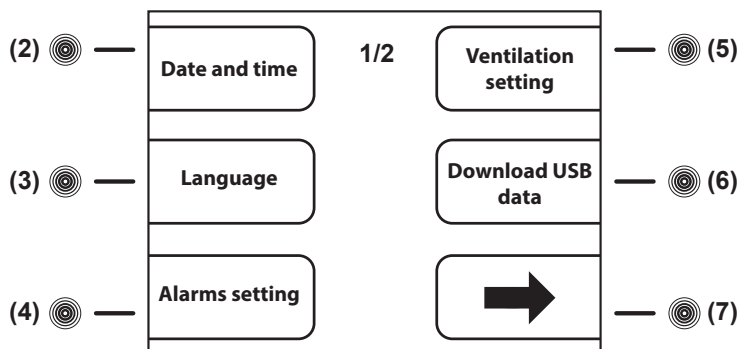
Fig.15

To access to the second page of the SETTING menu (**Fig.16**), press the **touch-button (7)→** from the first page of the SETTING menu (**Fig.15**).

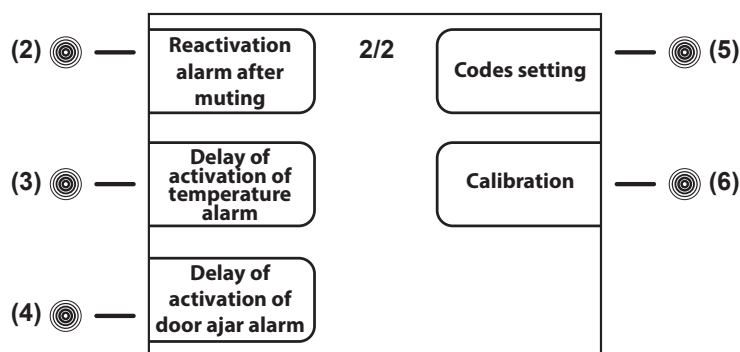
Fig.16

To set **Date and time** see paragraph 8.2.4.

8.3.6.1 Language setting (Fig.10)

To set the display language proceed as follow:


- deactivate the eventual touch-buttons locking protection (see par. 8.1.3)
- press the **MENU touch-button (9)** to access to the main menu (**Fig.12**)
- press the **touch-button (7)** to access to the setting menu (**Fig.15**)
- press the **touch-button (3)** to select the menu "Language"
- using the **touch-buttons (4)↑** and **(7)↓** to set the desired language and then press the **ENTER touch-button (12)** to confirm the new settings

8.3.6.2 Alarm setting (touch-button (4) Fig.15)

With this menu is possible to set the desired alarm limits for both the Chamber alarm limits (air temperature into the appliance) and the Product simulation alarm limits (temperature of the products stored into the appliance simulated with a ballasting of 250 ml of water or anti-freezing liquid).


To access to the ALARM SETTING menu, press the **touch-button (4)** while the display shows the first page of the setting menu (**Fig.15**).

Into the ALARM SETTING menu, there are present 2 sub-menus (**Fig.17**) that allow to modify the alarm limits both related to the chamber temperature and to the product simulation (if the optionally available extra sensor is installed).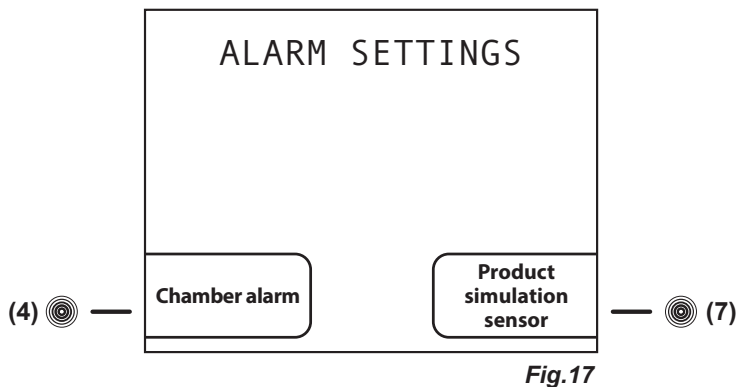

To set Chamber alarm limits, proceed as follow:


- from the sub-menu ALARM SETTINGS (**Fig.17**) press the **touch-button (4)** to access to the screen-shot CHAMBER ALARM SETTING (**Fig.18**).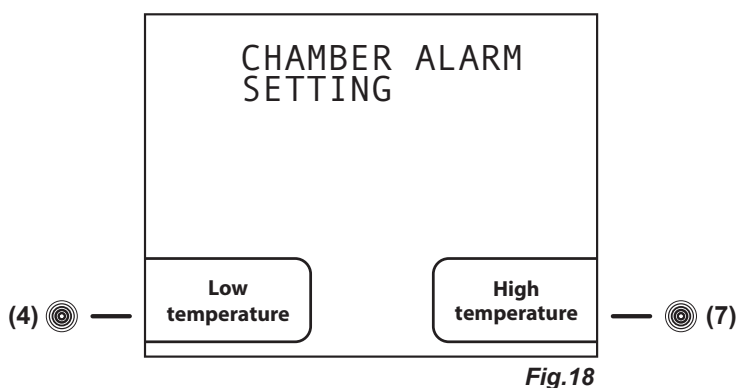

- press the **touch-button (4)** to access to the screen-shot SETTING CHAMBER LOW TEMPERATURE ALARM
- press the **touch-button (4)-** and **(7)+** to set the desired value
- or
- press the **touch-button (7)** to access to the screen-shot SETTING CHAMBER HIGH TEMPERATURE ALARM
- press the **touch-button (4)-** and **(7)+** to set the desired value
- press the **ENTER touch-button (12)** to confirm the new setting and the **ESC touch-button (11)** to return back to the previous menu or **HOME touch-button (10)** to return back to the main screen-shot

To set Product simulation sensor alarm limits, proceed as follow:

- from the sub-menu ALARM SETTING (**Fig.17**) press the **touch-button (7)** to access to the screen-shot PRODUCT ALARM SETTING (**Fig.19**)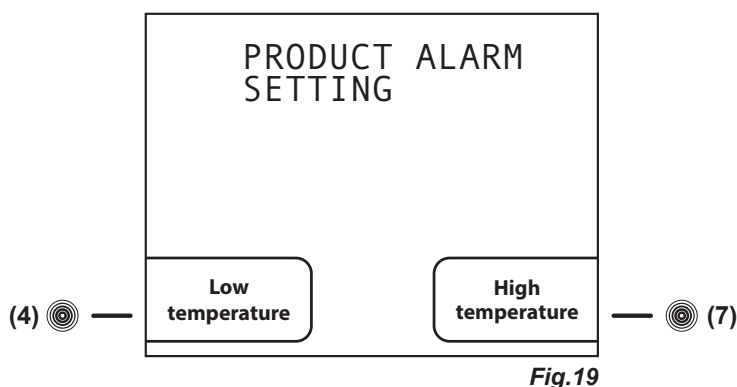

- press the **touch-button (4)** to access to the screen-shot SETTING PRODUCT LOW TEMPERATURE ALARM

- press the **touch-button (4)-** and **(7)+** to set the desired value
- or
- press the **touch-button (7)** to access to the screen-shot SETTING PRODUCT HIGH TEMPERATURE ALARM
- press the **touch-button (4)-** and **(7)+** to set the desired value
- press the **ENTER touch-button (12)** to confirm the new setting and the **ESC touch-button (11)** to return back to the previous menu or **HOME touch-button (10)** to return back to the main screen-shot

8.3.6.3 Ventilation setting (touch-button (5) Fig.15)

With this menu is possible to set the desired ventilation level into the appliance on 3 pre-fix levels.

To access to the VENTILATION SETTING, press the **touch-button (5)** while the display shows the first page of the setting menu (Fig.15).

- press the **touch-button (4)-** and **(7)+** to select among the 3 levels of ventilation available:
 - LOW: fan working in parallel with the compressor (apart door openings and defrost and some alarm conditions)
 - MEDIUM: fan working in parallel with the compressor and with spots functioning during the compressor OFF (apart door openings and defrost and some alarm conditions)
 - HIGH: fan working continuously (apart door openings and defrost and some alarm conditions)
- press the **ENTER touch-button (12)** to confirm the new setting and the **ESC touch-button (11)** to return back to the previous menu or **HOME touch-button (10)** to return back to the main screen-shot

8.3.6.4 Download USB data (touch-button (6) Fig.15)

With this menu (optionally available) is possible download on a USB memory-stick the data-logged details of the appliance, with the possibility to select the desired data to download. Following the instructions that will appear on the display, automatically will be downloaded a file (with .csv extension) on the USB memory-stick, that can be opened with MICROSOFT EXCEL[®] or whatever equivalent software compatible with the file extension. If present, the functionality, at the side of the xPRO control panel, there is a USB port.

The functionality can be even retro-fitted ordering the optionally available KIT USB EVERmed and become available after having correctly installed the software upgrade and the related accessories. The installation instructions are included in the KIT USB EVERmed. Once correctly installed the KIT USB EVERmed, it is possible to access to the Download USB data screen-shot and the subsequent functionalities.

To download the datas of interest, proceed as follow:

- To access to the DOWNLOAD USB DATA menu, press the **touch-button (6)** while the display shows the first page of the setting menu (Fig.15)
 - press the **touch-button (5)↑** and **(6)↓** to scroll among the list of the possible data to download
 - press the **touch-button (4)-** to clear the data (at the right of the display in correspondence of the data appears "no") or press the **touch-button (7)+** to select the data (at the right of the display in correspondence of the data appears "yes")
 - all the details with the indication "yes" are ready for the download
 - press the **ENTER touch-button (12)** to confirm and enter in the Download USB screen-shot
 - insert the USB memory-stick in the USB port
 - the display shows DOWNLOAD HISTORICAL DATA
 - press the **ENTER touch-button (12)** to confirm and enter in the DATE AND TIME
- Set the desired date and time from which download the datas. All the details will be automatically downloaded in the file. To set the date and time, proceed as follow:
- using the touch-buttons **(3)←** and **(6)→** highlight the position in the calendar or in the clock that should be modified
 - using the touch-buttons **(4)-** and **(7)+** to set the desired value
 - repeat the last 2 operations for all the numerical values in both calendar and clock, then press the **ENTER touch-button (12)** to confirm the details and process the download or the **ESC touch-button (11)** to return back to the previous menu
 - as soon will be finished the download, the display shows END-REMOVE USB
 - remove the USB memory-stick and check the file with a suitable PC

8.3.6.5 Reactivation alarm after muting (touch-button (2) Fig.16)

To access to the REACTIVATION ALARM AFTER MUTING, press the **touch-button (2)** while the display shows the second page of the setting menu (Fig.16).

- press the **touch-button (4)-** and **(7)+** to set the desired value (expressed in seconds, max. 900) that has to pass from the automatic reactivation of the acoustic alarm after the acoustic alarm signalling has been muted. The reactivation only works in case the alarm condition persists. If set the value "0", the reactivation is disabled
- press the **ENTER touch-button (12)** to confirm the new setting and the **ESC touch-button (11)** to return back to the previous menu or **HOME touch-button (10)** to return back to the main screen-shot

8.3.6.6 Temperature alarm activation delay (touch-button (3) Fig.16)

To access to the DELAY OF ACTIVATION OF TEMPERATURE ALARM, press the **touch-button (3)** while the display shows the second page of the setting menu (Fig.16).

- press the **touch-button (4)-** and **(7)+** to set the desired value (expressed in seconds, max. 900) that has to pass from the starting of the alarm condition of a temperature alarm till its signalling on the display and through the buzzer
- press the **ENTER touch-button (12)** to confirm the new setting and the **ESC touch-button (11)** to return back to the previous menu or **HOME touch-button (10)** to return back to the main screen-shot

8.3.6.7 Door ajar alarm activation delay (touch-button (4) Fig.16)


To access to the DELAY OF ACTIVATION OF DOOR AJAR ALARM, press the **touch-button (4)** while the display shows the second page of the setting menu (Fig.16).

- press the **touch-button (4)- and (7)+** to set the desired value (expressed In seconds, max. 600) that has to pass from the opening of the door till its signalling on the display and through the buzzer
- press the **ENTER touch-button (12)** to confirm the new setting and the **ESC touch-button (11)** to return back to the previous menu or **HOME touch-button (10)** to return back to the main screen-shot

8.3.6.8 Codes setting (touch-button (5) Fig.16)

From this menu is possible to customize the 2 appliance access codes. The **ON/OFF code**, which is required every time that it is needed to switch the appliance ON or OFF (**factory code 0050**). The **Keyboard blocking code**, which has to be entered every time that the appliance has been unused for 5 minutes (**factory code 0010**).

To access to the screen-shot ACCESS CODE SETTING (Fig.20), press the **touch-button (5)** while the display shows the second page of the setting menu (Fig.16).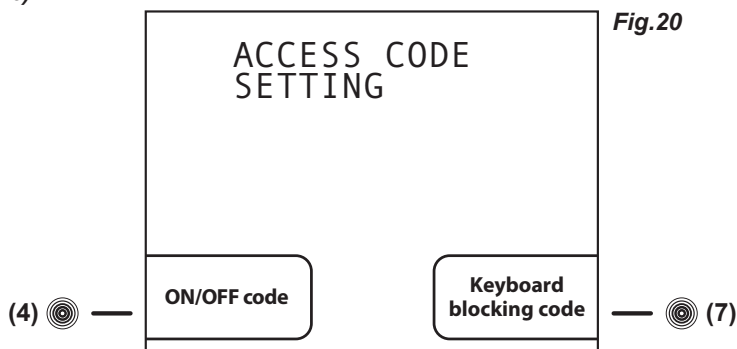

To set a new ON/OFF code, proceed as follow:

- press the **touch-button (4)** to access to the screen-shot ON/OFF CODE
- using the touch-buttons **(3)←** and **(6)→** highlight the position in the code that should be modified
- using the touch-buttons **(4)-** and **(7)+** to set the desired value
- press the **ENTER touch-button (12)** to confirm the new setting and the **ESC touch-button (11)** to return back to the previous menu


To set a new Keyboard blocking code, proceed as follow:

- press the **touch-button (4)** to access to the screen-shot KEYBOARD BLOCKING CODE
- using the touch-buttons **(3)←** and **(6)→** highlight the position in the code that should be modified
- using the touch-buttons **(4)-** and **(7)+** to set the desired value
- press the **ENTER touch-button (12)** to confirm the new setting and the **ESC touch-button (11)** to return back to the previous menu

NB: Setting the value "9999" function will be disabled.

8.3.6.9 Calibration (touch-button (6) Fig.16)

From this menu is possible to manually calibrate 2 sensors of the appliance. The Chamber sensor is the one acquiring the temperature into the appliance, while the Product simulation sensor (optionally available) simulate the temperature of the products stored into the appliance with a ballasting of 250 ml of water (or anti-freezing liquid). Through this menu is possible to adjust the reading of the sensors according to reading of external comparison thermometers, data loggers, thermocouples or personal wishes. To access to the CALIBRATION SETTING (Fig.21), press the **touch-button (6)** while the display shows the second page of the setting menu (Fig.16)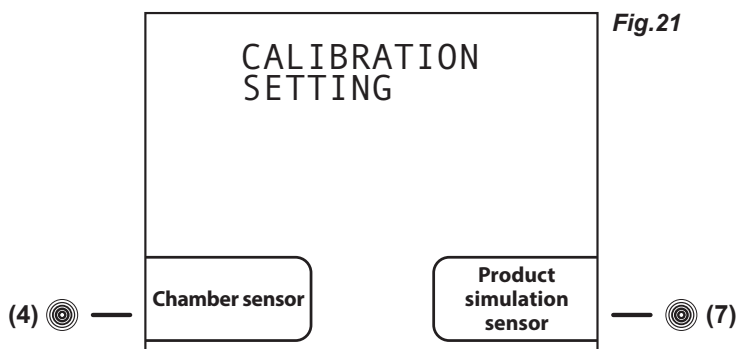

To set a new calibration for the Chamber sensor, proceed as follow:

- press the **touch-button (4)** to access to the screen-shot CALIBRATION CHAMBER SENSOR
- press the **touch-button (4)- and (7)+ to increase** (max + 10°C) or lower (min -10°C) the calibration value highlighted
- press the **ENTER touch-button (12)** to confirm the new setting and the **ESC touch-button (11)** to return back to the previous menu or **HOME touch-button (10)** to return back to the main screen-shot

To set a new calibration for the Product simulation sensor (optionally available), proceed as follow:

- press the **touch-button (4)** to access to the screen-shot CALIBRATION PRODUCT SENSOR
- press the **touch-button (4)- and (7)+ to increase** (max +10°C) or lower (min -10°C) the calibration value highlighted
- press the **ENTER touch-button (12)** to confirm the new setting and the **ESC touch-button (11)** to return back to the previous menu or **HOME touch-button (10)** to return back to the main screen-shot

8.4 OPERATION OF THE TEMPERATURE CHART RECORDER


8.4.1 Description of the temperature chart recorder (Fig.22)

The appliance may be fitted (present if ordered in some series), with the temperature chart recorder, recording on paper charts the appliance internal temperature. The temperature recorder is available in different temperature ranges according to the model on which it is installed:

- Range : -10°C / +40°C
- Range : -35°C / +15°C
- Range : -50°C / +50°C


The survey of the chamber temperature is done with a sensor independent from the control panel sensor, while the recording, weekly operation, is on paper charts with an ink-tip. The temperature recording is granted even during power failure periods thanks to the battery powering of the recording device. The battery is 1,5V, AA type. A perspex cover allows checking the diagrams but at the same time it protects them against tampering of unauthorized personnel thanks to the key locks safety.

Fig.22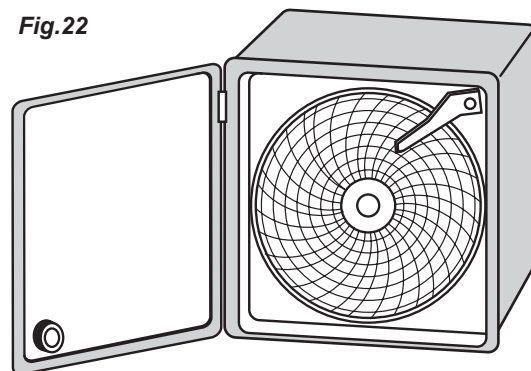


8.4.1.1 Paper chart replacement

- Open the perspex cover
- Unscrew the fixing nut
- Lift gently the metal arm on which is fitted the ink-tip
- Remove the installed paper chart, paying attention at the 2 metal edges surrounding the chart that keep it in position
- Insert the new paper chart, paying attention to insert it into the central hinge and into the 2 metal edges surrounding the chart
- Lower gently the metal arm on which is fitted the ink-tip in order to return it in the original position
- Drive the paper chart in order to position the ink-tip writing point in the exact point from which it has to start to record the temperature, paying attention to do not write on the diagram. In order to choose the exact recording start point, refer to the days and times printed on the chart itself
- Screw the fixing nut
- Close the perspex cover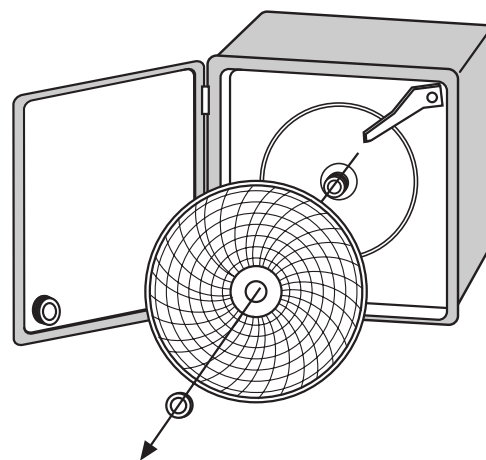


8.4.1.2 Ink-tip replacement

- Open the perspex cover
- Lift gently the metal arm on which is fitted the ink-tip
- Remove, pulling, the ink-tip from the metal arm
- Insert the new ink-tip on the metal arm paying attention to the guideways on the ink-tip itself. Push until reach the limit-stop
- Lower gently the metal arm on which is fitted the ink-tip in order to return it in the original position
- Close the perspex cover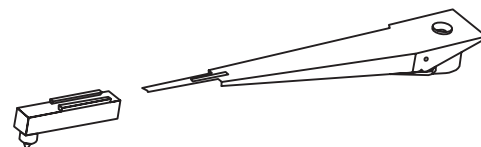

8.4.1.3 Battery replacement

- Open the perspex cover
- Unscrew the fixing nut
- Lift gently the metal arm on which is fitted the ink-tip
- Remove the installed paper chart, paying attention at the 2 metal edges surrounding the chart that keep it in position
- Pull the central hinge on which was screwed the nut, and at the same time moving slightly up and down to make easier the extraction of the clockwise mechanism
- Replace the battery in the rear side of the clockwise mechanism paying attention to the battery polarity
- Insert the clockwise mechanism in the hole left on the chart recorder, simply making pressure on the mechanism
- Insert the paper chart, paying attention to insert it into the central hinge and into the 2 metal edges surrounding the chart
- Lower gently the metal arm on which is fitted the ink-tip in order to return it in the original position
- Drive the paper chart in order to position the ink-tip writing point in the exact point from which it has to start to record the temperature, paying attention to do not write on the diagram. In order to choose the exact recording start point, refer to the days and times printed on the chart itself
- Screw the fixing nut
- Close the perspex cover

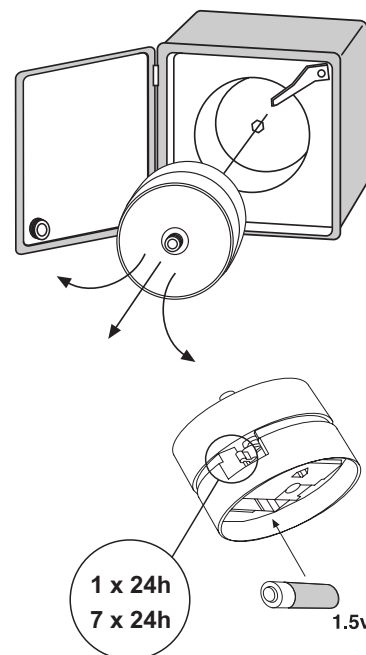8.5 OPERATION OF THE DIGITAL PRINTER

8.5.1 Description of the digital printer

The appliance may be fitted (present if ordered in some series), with the temperature digital printer.

The digital printer records the temperature in an internal non-volatile memory the temperature of the appliance. The recording periods and intervals are fully adjustable and can be easily set through the printer soft-touchpads and the wide display (with clear written indications). Then, the temperature trends can be printed on paper, selecting among 3 printing modes according to the need, online (printing out at fix intervals), on daily basis (once per day) or an historical report of all the datas acquired.

The survey of the chamber temperature is done with a sensor independent from the control panel sensor. To increase the safety level, it is included an alarm for sensor failure. The printing is ink-free on thermal paper.

For more detailed information, please refer to the specific booklet (included among the instructions if the printer is present).

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EVERmed Srl: Via Galileo Galilei, n° 2 - 46020 Motteggiana (MN) ITALY